

Product bulletin

ChemSeal™

Alfa Aesar's ChemSeal™ packaging is specifically designed for air, oxygen and moisture sensitive and volatile compounds. ChemSeal packaging allows the user to sample liquids through the resealable septum. Using a syringe, a volumetric quantity of the reagent can be sampled and transferred to your reaction medium without exposure to moisture or oxygen. Upon withdrawing the syringe, ChemSeal reseals the bottle. The end result is a better conservation of your stock reagent and a better control over your chemicals. Additional seals are also available.

ChemSeal usage instructions

Please follow these instructions for safe and efficient use of our ChemSeal product:

- Do not unscrew the cap.
- Use appropriately sized syringe to withdraw the required sample.
- Keep the product under argon.

